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INSTRUCTION MANUAL

ZOOM STEREO MICROSCOPE XTS-SP2

Thank you for purchasing the Zoom Stereo Microscope.
Please read this instruction manual thoroughly before use to insure proper operation and long service.


NOTES

- This is a precision instrument – Handle With Care. Use caution to avoid shock or excessive forces on instrument when in use or transport.
- Avoid direct sunlight or areas with high temperatures or high levels of dust. Keep in location where temperate and humidity are stable.
- Keep lens surface clean from dirt and fingerprints.
- Clean the lens gently with lens cleaning cloth or soft cotton cloth.
- Do not turn left and right Zoom Controls in opposite directions – doing so will result in damage to instrument.
- When not in use, store with protective cover, in location that has low moisture in order to prevent internal mold growth.

SPECIFICATIONS

Magnification	6.5~45×・13~90×
Objective Lens	0.65~4.5× (Zoom Ratio 1:7)
Eyepiece Lens	WF10× (20mm Field of View) WF20× (10mm Field of View)
Working Distance	95mm
Eyepiece Angle	45°
Focus Adjustment	Movable Body
Stand	Pole Type
Stage Plate	φ 95 Black/White Acrylic
Eye Width Adj. Range	55~75mm
Diopter Adj. Range	± 5 Diopter
Illumination	Optional Ring Light (sold separately)
Ring Light Attachment Dimensions	Light Mounting Section : φ 54 mm Microscope Mounting Threads : M48×P0.75
Base Dimensions	L245×W170×H40mm
Weight	Approx. 5.4 kg

PART IDENTIFICATION - ASSEMBLY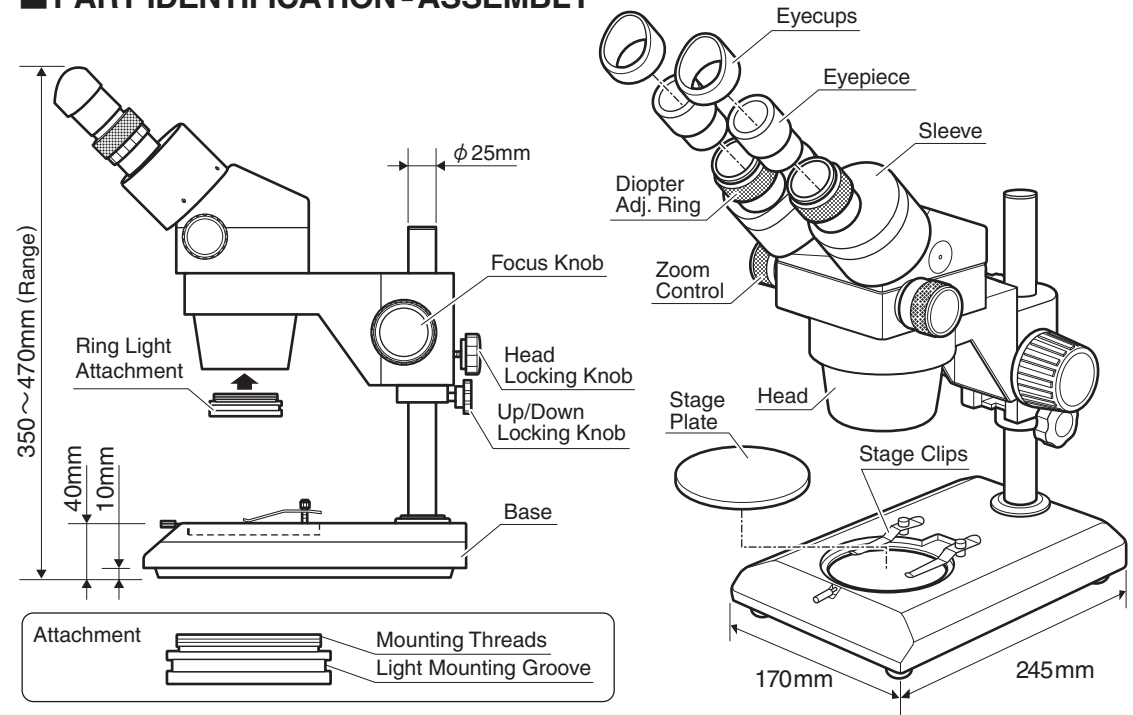

OPERATION

1. Sample Placement

- ① Place sample in center of Stage.
- ② If required, use Stage Clips to hold sample in place.

2. Eyecup Use

- **For use without eyeglasses:**
Hold Diopter Adjustment Ring and rotate Eyepiece to adjust diopter setting. Align the Eyecups to point toward user and look through the Eyepiece with eyes brought close.
- **For use with eyeglasses:**
Remove Eyecups to use.

3. Eye Width Adjustment

Hold the left and right sleeve in each hand and look through the Eyepieces. Adjust the Sleeves to align the left and right field of view seen through the Eyepieces.

4. Focus Adjustment • Diopter Adjustment

- ① Rotate the Zoom Control for highest magnification.
- ② Look through the right eyepiece with the right eye and adjust the focus using the Focus Knob.
- ③ Rotate the Zoom Knob for lowest magnification.
- ④ Using the Right Diopter adjustment, adjust the focus if required.
- ⑤ Now rotate the Zoom Control back to highest magnification and check the focus. Repeat steps 2 to 4 to achieve the highest focus accuracy.
- ⑥ Rotate the Zoom Control for highest magnification. Look through the left Eyepiece with the left eye, and adjust focus using the Left Diopter Adjustment Ring.

SC Niigata seiki Co., Ltd.

5-3-14, Tsukanome, Sanjo, Niigata, Japan, 955-0055
Tel. : +81-256-33-5522 Fax. : +81-256-33-5518
MAIL intl.sales@niigataseiki.co.jp
URL http://www.niigataseiki.co.jp