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

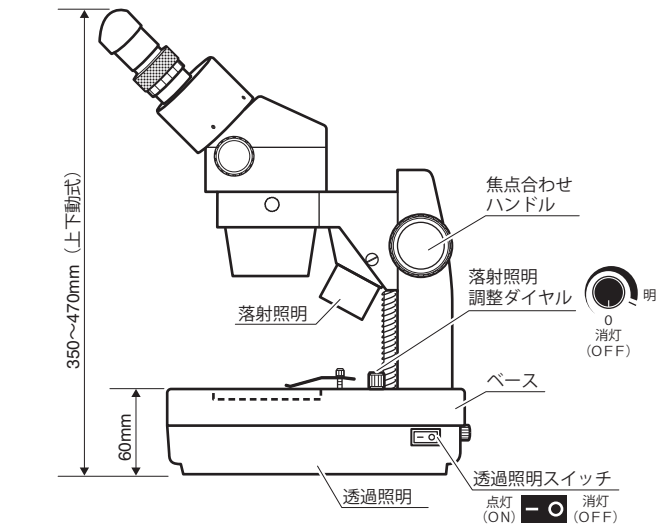
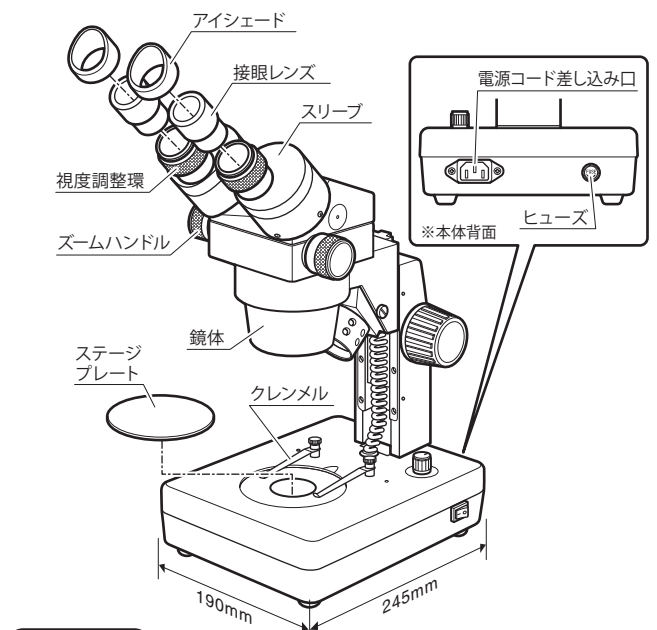
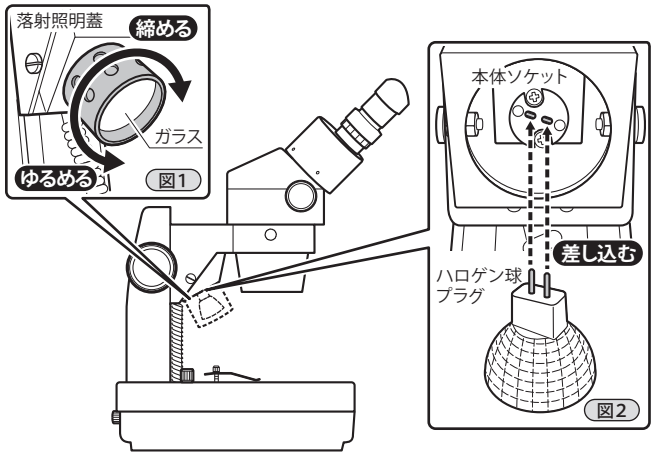
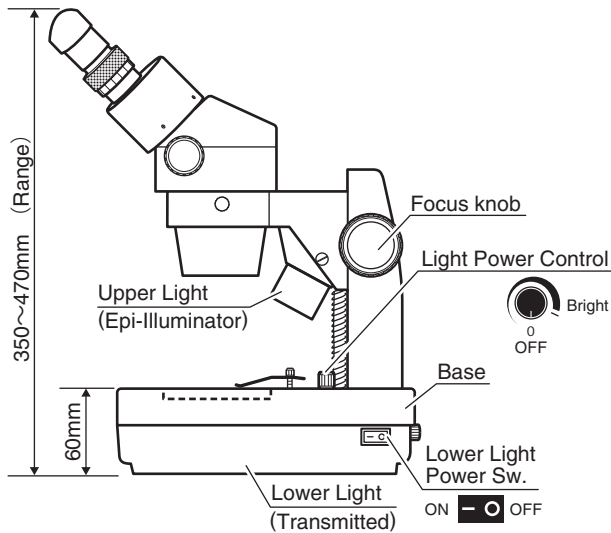
STEREO ZOOM MICROSCOPE

Thank you for purchasing the Stereo Zoom Microscope.
Please read this instruction manual thoroughly before use to insure proper operation and long service.

PART IDENTIFICATION & SETUP

<Bulb Installation>

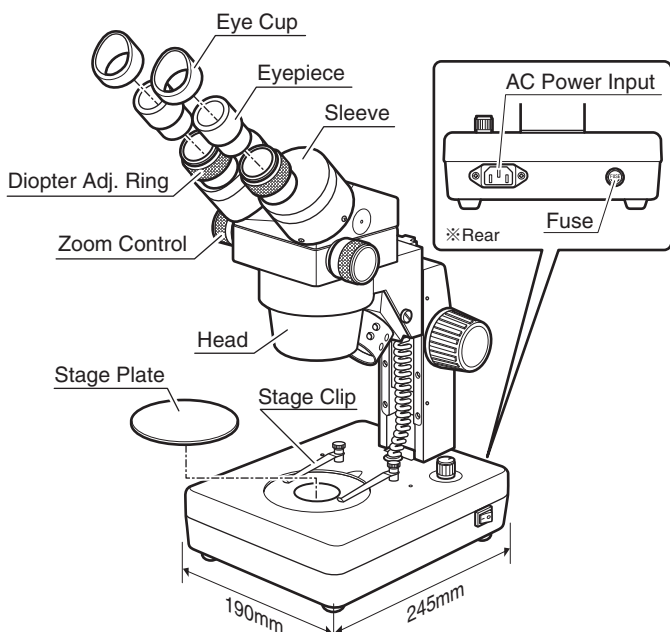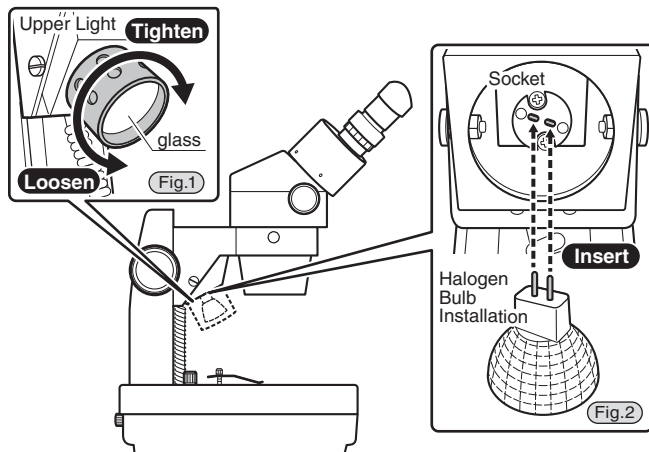⚠ CAUTION

- Only use as directed. Improper use may cause accident or injury.
- When not in use store with protective cover, in location that has low moisture in order to prevent internal mold growth.
- Lighting unit not certified for use outside of Japan.

SPECIFICATIONS

Rated Voltage	AC100V 50/60Hz
Magnification	6.5~45× (for 10× eyepiece)
	13~90× (for 20× eyepiece)
Objective Lens	0.65~4.5× (Zoom Ratio 1:7)
Eyepiece Lens	WF10× (FOV = 20mm) · WF20× (FOV = 10mm)
Working Distance	95mm
Eyepiece Angle	45°
Focus Adjustment	Movable Body
Stand	Rack and Pinion Type
Stage Plate	φ95mm×2pcs ①Ground Glass ②black/white (plastic)
Eye Width Adj. Range	55~75mm
Diopter Adj. Range	±5diopter
Illumination	Epi-illuminator: 6V·15W Halogen bulb
	Transmitted: 5W ring light
Base	L245×W190×H60mm
Weight	approx.5.7kg

NOTES

- This is a precision instrument – Handle With Care. Use caution to avoid shock or excessive forces on instrument when in use or transport.
- Use and store in location with low humidity and stable temperatures. Avoid locations with dust, direct sunlight, excessive heat, or high vibration.
- Keep lens surface clean from dirt and fingerprints.
- Clean the lens gently with lens cloth or soft cotton cloth.
- Do not turn the left and right Zoom Controls in opposite directions; this will damage the instrument.
- Before use, install the halogen bulb following the procedure below.

HALOGEN BULB INSTALLATION

- ① Remove the bulb cover from the Upper Light by rotating as shown. Use care not to drop or damage glass. (Fig.1)
- ② Insert the bulb into the 2-prong socket. (Fig.2)
 - ※ If not fully inserted, the bulb will not light but use care not to force or damage bulb when inserting.
- ③ Replace the bulb cover and rotate to tighten. (Fig.1)

SAMPLE PLACEMENT

- ① Place the sample in the center of the Stage Plate.
- ② If required, use Stage Clips to hold sample in place.

EYE CUP USE

- For use without eyeglasses: Hold Eyepiece and rotate the Diopter Adjustment Ring to adjust diopter setting. Align the Eye cups to point toward user and look through the Eyepiece with eyes brought close.
- For use with eyeglasses: Remove Eye cups to use.

EYE WIDTH ADJUSTMENT

Hold both sleeves in each hand and look through the Eyepieces. Adjust the Sleeves to align the left and right field of view seen through the Eyepieces.

FOCUS & DIOPTER ADJUSTMENT

- ① Rotate the Zoom Control for highest magnification.
- ② Look through the right eyepiece with the right eye and adjust the focus using the Focus Knob.
- ③ Rotate the Zoom Knob for lowest magnification.
- ④ Using the Right Diopter adjustment, adjust the focus if required.
- ⑤ Now rotate the Zoom Control back to highest magnification and check the focus. Repeat steps ②~④ to achieve the best focus accuracy.
- ⑥ Rotate the Zoom Control for highest magnification. Look through the left Eyepiece with the left eye, and adjust focus using the Left Diopter Adjustment Ring.

FUSE REPLACEMENT

In case of fuse failure, it can be replaced with a commercially available 250V 2A fuse.