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MEMO

Thank you for purchasing the USB Microscope Camera. Please read this instruction manual thoroughly before use to insure proper operation and long service.

FEATURES

- 1.3M pixel USB Camera designed for easy use on a microscope. Simply connect to USB port for video image display.
- Powered by USB port; No separate power supply required.

NOTES


- This is a precision instrument - Handle With Care. Use caution to avoid shock or excessive forces on instrument when in use or transport.
- Do not use or store in direct sunlight or areas with high temperatures or high levels of dust. Keep in location where temperature and humidity are stable.
- Keep lens surface clean from dirt and fingerprints. To clean the lens, wipe gently with lens cleaning cloth or soft cotton cloth.
- When not in use, store in supplied case in location which has low moisture in order to prevent corrosion or internal mold growth.

SPECIFICATIONS

Model	MU-130
Sensor	1/3.2 CMOS
Pixel Count	130M Pixels
Frame Rate	30Frames/sec.
Interface	USB 2.0
Power Supply	Supplied by USB port
Supported OS	Windows XP (SP3), Vista, 7, 8, 10
Min. recommended PC requirements	<ul style="list-style-type: none"> ● CPU: 2.4GHz clock ● Memory: 1GB ● Hard Drive: 10GB Free space
Weight (Body + Lens)	approx.175g
Accessories	<ul style="list-style-type: none"> ● Software CD ... 1 ● Bore Adapters: <ul style="list-style-type: none"> φ 23.2→φ 30.0mm ... 1 φ 23.2→φ 30.5mm ... 1 ● USB Cable (-1.5m) ... 1

1 PART NAMES & ASSEMBLY

- 1 Remove the eyepiece from the microscope.
- 2 Remove the Lens Caps from the Camera Body and the Lens.
- 3 Follow the diagram below to connect.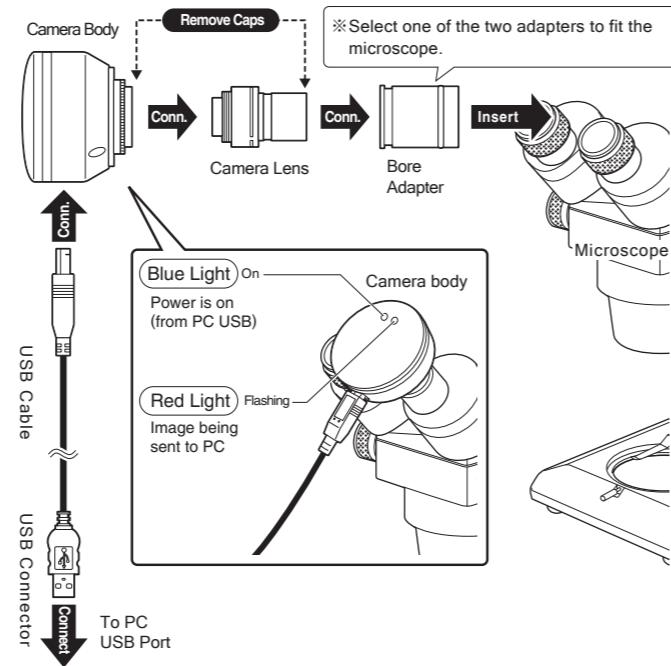


- 4 When MU-130 Camera is connected to the PC USB port using the cable, the following message will appear in the bottom right corner of screen. Please proceed.
- ※ Screen appearance will vary for different PC settings.
 ※ Keep the MU-130 camera connected to the USB port and continue with the additional installation steps. Keep the MU-130 camera connected to the USB port and continue with the additional installation steps.


2 INSTALLING ScopelImage 9.0 SOFTWARE

Administrator privileges for computer are required for installation.


- 1 Double click on the file "SISetupX1N.exe" in the CD-ROM folder.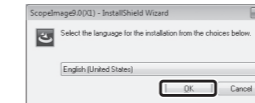


- 2 When the language selection screen appears, select "English" and click "OK"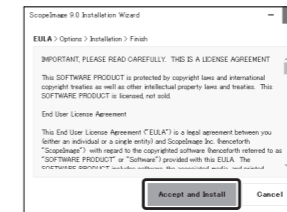

- 3 When the installation guide screen appears, click on "Accept and Install" to continue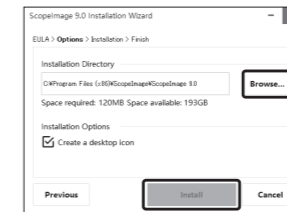

- 4 Click on the "Browse" button to specify the directory for installation; when done click on "Install" to start the installation.

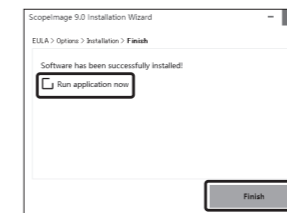- 5 Click on "Finish" when the installation is complete.
 ※ To start the software immediately once installation is complete, check the box labeled "Run application now"

- 6 The icon shown to the right will be placed on the desktop when the viewer software is installed.

3 USING THE VIEWER SOFTWARE (ScopelImage 9.0)


Feature Support

Viewer Software (ScopelImage 9.0) provides supports for the following features:

- Image preview display
 - Preview screen control toolbar
 - Still image capture and save
 - Video capture and save
- Other features are not covered. Please note, support for additional features can not be provided even if requested.

1 Image Preview


- Double click on ScopelImage9 icon to start the application.
- ※ Select the device labeled "Camera-X1N"
- Using the preview screen, follow the microscope instructions for adjusting focus of the object being viewed.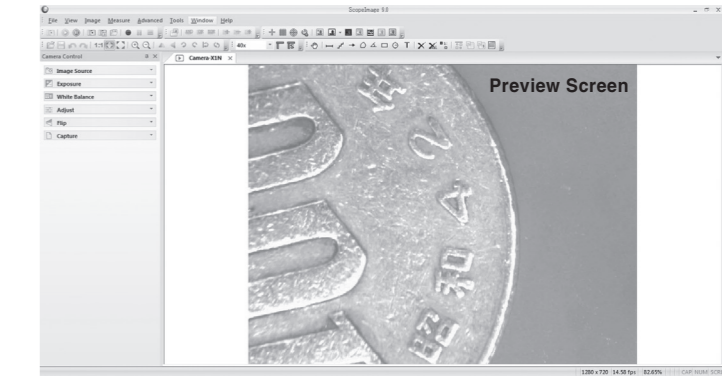

3 USING THE VIEWER SOFTWARE (ScopelImage 9.0) - continued

2 Image Preview Operation

- The function toolbar can be used to manipulate the image displayed on the preview screen.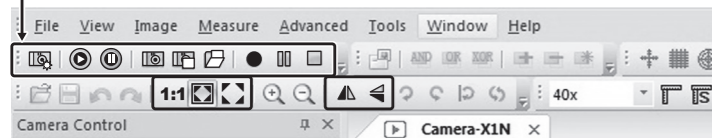


	Open preview screen. When icon is not lit, preview can not be opened.	1:1	Display pixels in without scaling for highest details when adjusting focus.
	Play or resume the preview screen.		Display full field of view for adjusting size of image shown on preview screen.
	Pause the preview screen display.		Full screen display on monitor (click on "Close Full Screen" button to close).
	Save the displayed preview as a still image. (→[3])		Flip the preview display along vertical axis (right and left.) ※ 1
	Freeze the image displayed on the preview screen. ※ To save image, click on: File menu → Save Clicking on the [X] in the image tab closes the image.		Flip the preview display along horizontal axis (up and down.) ※ 1
	Open the destination folder.		
	Save a video of the image. (→[4])		
	Pause video capture. (→[4])		
	Resume video capture. (→[4])		

※1 Depending on PC settings, these features may not be operable.


3 Still Image Capture and Save

- The preview screen image can be frozen and saved as an image file by clicking on the image capture icon on the toolbar.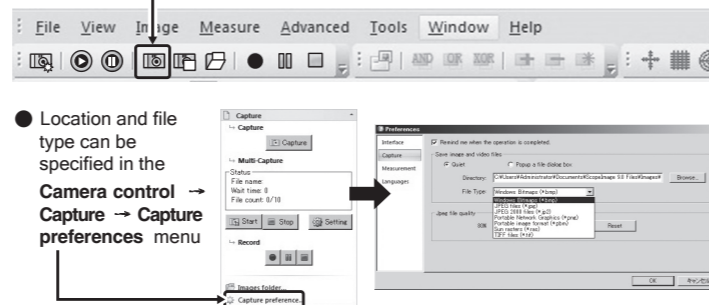


4 Video Capture and Save

- Image displayed on the Preview Screen can be captured and saved as video.

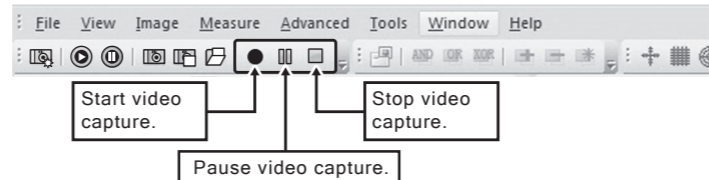- Image format can be specified in the **Camera control → Image Source** menu.

- Select the image format based on your requirements:
- ◆ MJPG Select when high frame rate is desired.
 - ◆ YUY2 Select for high image quality.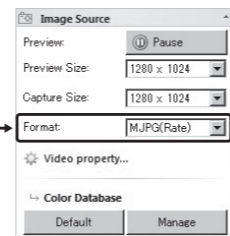

4 UNINSTALL PROCEDURE ※ Will vary for different PC settings

- In the Windows Control Panel, Select "Uninstall a program" and then select ScopelImage9.0 and uninstall.

5 TROUBLESHOOTING

- Q. Display image is dark or blurry.**
 - A. Insure there is sufficient lighting using microscope lamp or accessory light.
- Q. Driver software does not install. Viewer software does not work properly.**
 - A. Please confirm PC operating system is compatible; contact distributor.
- Q. Shortcut does not appear on the desktop.**
 - A. Shortcut will only be created for user with administrator privileges. If you want to create a shortcut for user without administrator privileges, go to the application installation directory to create the shortcut.

Note) For default installation, the application is installed in the following directories: C:\Program Files\ScopeImage\ScopImage 9.0
 C:\Program Files (x86)\ScopeImage\ScopImage 9.0

MEMO