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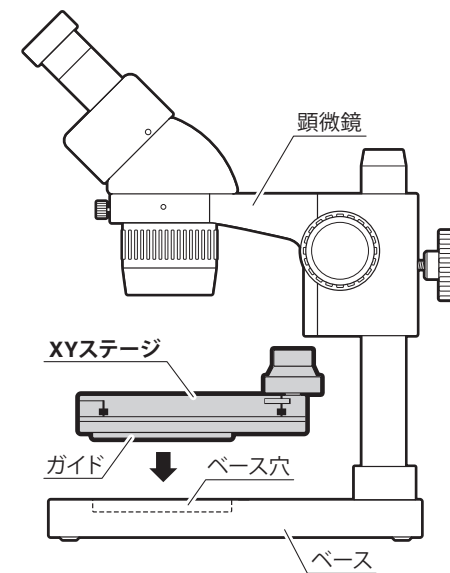
INSTRUCTION MANUAL

XY MICROSCOPE STAGE STD-8090

Thank you for purchasing Microscope X-Y STAGE STD-8090.
Please read this instruction manual thoroughly before use to insure proper operation and long service.

APPLICATION

- Placed upon base of microscope, fine adjustment of sample horizontal and vertical position can be smoothly controlled using the adjustment dial.

NOTES


- This is a precision instrument. Avoid excessive shock or pressure when transporting and during use. Please handle with care.
- Avoid using or storing in locations where there is a lot of dust, direct sunlight, or subject to high temperatures. Please chose a location with low humidity and temperature changes.
- Glass stage can scratch easily. Wipe lightly using lens cleaning cloth only to clean.
- Water or dust entering the internal controls can cause failure. Please handle sample carefully to avoid damaging the stage.
- Do not disassemble unit.
Do not remove screws other than removable guide mounting screws.
- Store in low humidity conditions to avoid rust and mold.

SPECIFICATIONS


Model No.	STD - 8090
Maximum sample load	5.5 kg
Dimensions	approx. 180 x 175 x 57mm (including adjustment dial)
Glass stage dimensions	105 x 143mm (Vertical x Horizontal)
Stage range of motion	80 x 90mm (Vertical x Horizontal)
Compatible microscope base	φ95mm
Material	Body : Aluminum, Stage : Glass Adjustment dial : Plastic
Weight	approx. 1.7 kg
Accessories	No-slip sheet (rubber) x 1

SET UP AND PARTS IDENTIFICATION

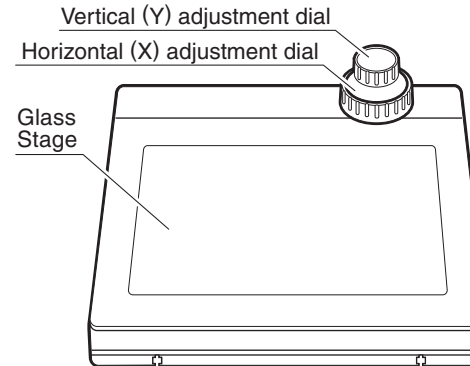TOP

BOTTOM with No-slip sheet

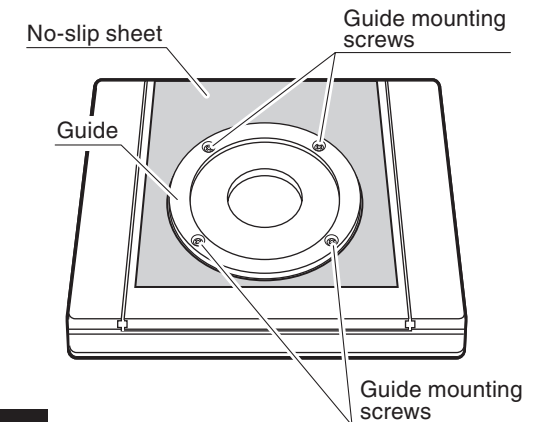
1. No-slip sheet installation

※If No-slip sheet is required, please use pressure sensitive adhesive to apply to the bottom of the stage.

- Peel off the white release paper from No-slip sheet.
- Apply No-slip sheet to bottom of stage as shown in the illustration.

2. Installation of the X-Y Stage onto microscope

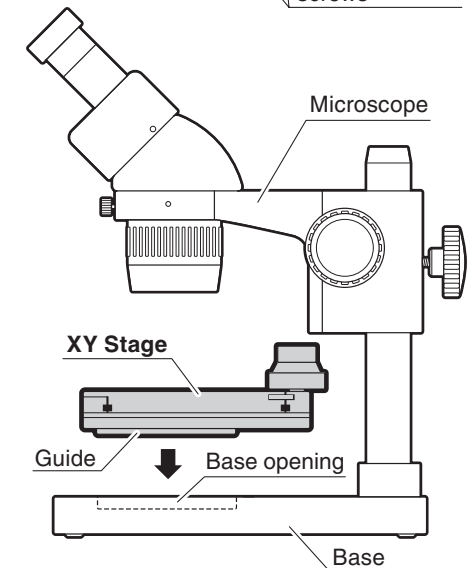- Remove specimen clips and sample plate from base of microscope.
- Using the guide on the bottom of the stage, set the stage into the opening in the base of the microscope.

For microscopes with non-compatible base openings

※The guide on the X-Y stage can be removed for use when the stage does not fit the diameter of the microscope base opening.

- Remove the guide mounting screws using a #2 Phillips driver (4 locations.)
- Place the X-Y Stage on the microscope base.

OPERATION

- Place the specimen on the stage glass.
- While looking through the microscope, use the adjustment dial to fine tune the sample position.

SK Niigata Seiki Co., Ltd.

5-3-14, Tsukanome, Sanjo, Niigata, Japan, 955-0055
Tel. : +81-256-33-5522 Fax. : +81-256-33-5518
MAIL intl.sales@niigataseiki.co.jp
URL <https://www.niigataseiki.net/en/>