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## INSTRUCTION MANUAL

# ROTARY STAGE FOR MICROSCOPE STD-360XY

Thank you for purchasing Rotary Stage For Microscope.  
Please read this instruction manual thoroughly before use to insure proper operation and long service.

### APPLICATION

- Place upon Microscope Base for precise control of sample position. Adjustment Dials adjust sample position Right-Left, Front-Back, and Rotation.

### NOTES


- This is a precision instrument. Avoid excessive shock or force when transporting and during use. Please handle with care.
- Avoid using or storing in locations where there is a lot of dust, direct sunlight, or subject to high temperatures. Please chose a location with low humidity and temperature changes.
- Glass stage can scratch easily. Wipe lightly using lens cleaning cloth only to clean.
- Water or dust entering the internal controls can cause failure. Please handle sample carefully to avoid damaging the stage.
- Do not disassemble unit.  
Do not remove screws other than removable guide mounting screws.
- Store in low humidity conditions to avoid rust and mold.

### SPECIFICATIONS


Model No.	STD-360XY
Maximum sample load	4.0kg
Dimensions	approx. 180×165×65mm (including adjustment dial)
Glass stage dimensions	approx. $\phi$ 95mm
Stage range of motion	Y:~55mm • X:~75mm • Rotation:360°
Compatible microscope base	approx. $\phi$ 95mm
Material	Body : Aluminum, Stage : Glass Adjustment dial : Plastic • Aluminum
Weight	approx. 2kg
Accessories	No-slip sheet (rubber) x 1

### SET UP AND PARTS IDENTIFICATION

【TOP】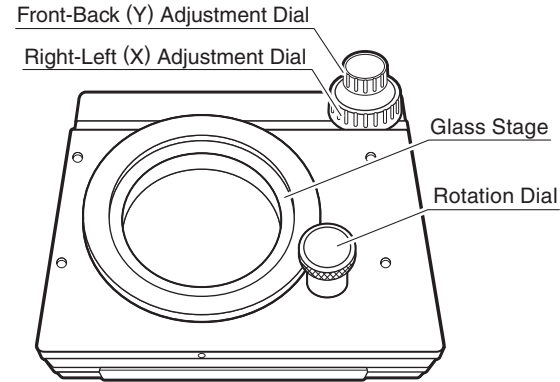

【BOTTOM】 with No-slip sheet

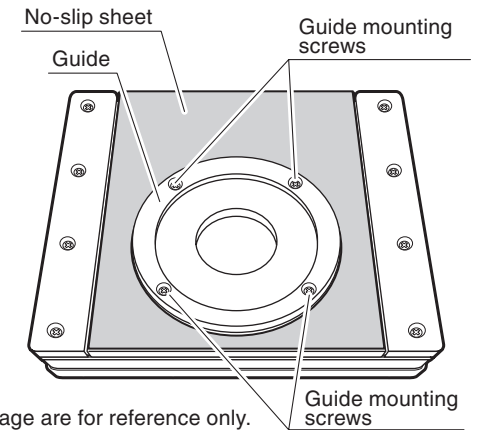※The distance scales on the Adjustment Dials and X-Y Stage are for reference only.

### Installation of the X-Y Stage onto microscope

- ① Remove specimen clips and sample plate from base of microscope.
- ② Using the guide on the bottom of the stage, set the stage into the opening in the base of the microscope.
- ③ Place the Glass Stage onto the main stage.

**CAUTION** The Glass Stage is not attached to the main stage. Use care not to drop when setting up or transporting.

### No-Slip Sheet


※If No-slip sheet is required, please use pressure sensitive adhesive to apply to the bottom of the X-Y stage.

- ① Peel off the white release paper from No-slip sheet.
- ② Apply No-slip sheet to bottom of X-Y stage as shown in the illustration.

### For microscopes with non-compatible base openings

※If base opening is too shallow, applying the No-slip sheet may improve the fit.  
Also, if the Guide causes interference, it can be removed by following these steps.

- ① Remove the guide mounting screws on the underside of the stage using a #2 Phillips driver (4 locations.)
- ② Place the X-Y Stage on the microscope base.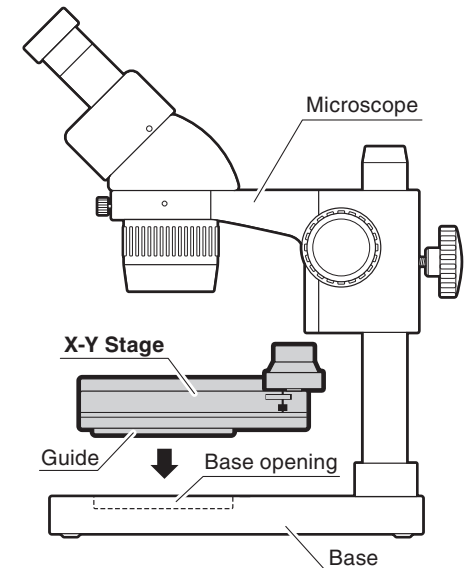

### OPERATION

- ① Place the specimen on the stage glass.
- ② While looking through the microscope, use the adjustment dials to fine tune the sample position.

**Niigata Seiki Co., Ltd.**

5-3-14, Tsukanome, Sanjo, Niigata, Japan, 955-0055  
Tel. : +81-256-33-5522 Fax. : +81-256-33-5518  
MAIL intl.sales@niigataseiki.co.jp  
URL https://www.niigataseiki.co.jp