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INSTRUCTION MANUAL

ZOOM STEREO MICROSCOPE XZ-45N

Thank you for purchasing the NSK Zoom Stereo Microscope. Please read this instruction manual thoroughly before use to insure proper operation and long service.


■NOTES

- This is a precision instrument – Handle With Care. Use caution to avoid shock or excessive forces on instrument when in use or transport.
- Avoid direct sunlight or areas with high temperatures or high levels of dust. Keep in location where temperate and humidity are stable.
- Keep lens surface clean from dirt and fingerprints.
- Clean the lens gently with lens cleaning cloth or soft cotton cloth.
- Do not turn the left and right Zoom Controls in opposite directions; this will damage the instrument.
- When not in use store with protective cover, in location that has low moisture in order to prevent internal mold growth.

■SPECIFICATIONS

Magnification	10～45×
Objective Lens	1～4.5× (Zoom Ratio 1:4.5)
Eyepiece Lens	WF 10× (20mm Field of View)
Working Distance	97mm
Eyepiece Angle	60°
Focus Adjustment	Movable Body
Stand	Pole Type
Stage Plate	φ95mm Black/White Acrylic
Diopter Adj. Range	±5Diopter
Illumination	Optional Ring Light (sold separately) Mounting diameter: φ53mm
Base Dimensions	L 255×W 200×H 25mm
Weight	approx. 3.6kg
Accessories	・Eyecup...2 ・Protective Cover...1

■PART IDENTIFICATION and ASSEMBLY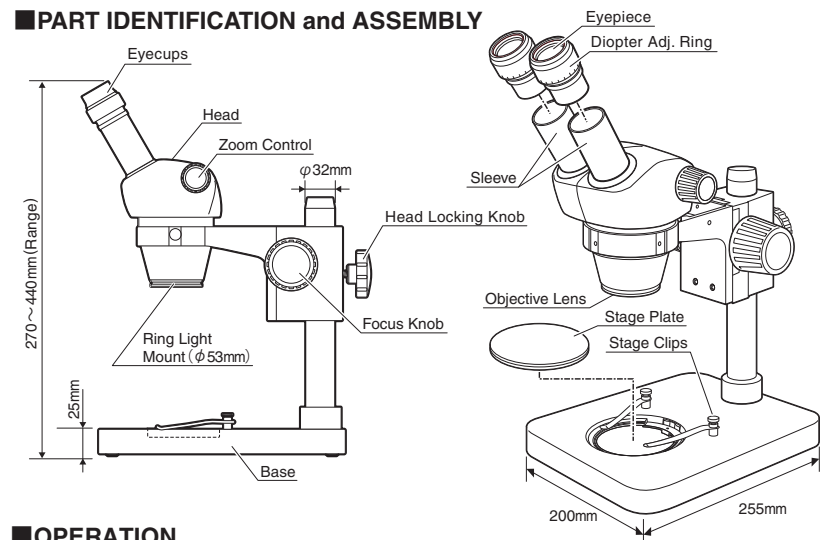

■OPERATION

1. Sample Placement


- ① Place the sample in the center of the Stage Plate.
- ② If required, use Stage Clips to hold sample in place.

2. Eyecup Use

- ① Remove the rubber Guard from the Eyepieces.
- ② Install the Eyecups as shown.

※ Make sure the Eyecups are seated into the groove on the Eyepiece.

※ When wearing eyeglasses, keep the Guards in place and do not install the Eyecups.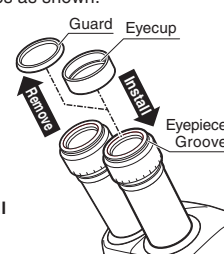

3. Focus Adjustment • Diopter Adjustment

- ① Rotate the Zoom Control to set for highest magnification.
- ② With right eye, look through the right eyepiece and adjust the focus using the Focus Knob.
- ③ Rotate the Zoom Control for lowest magnification.
- ④ Using the Right Diopter Adj. Ring, adjust the focus if required.
- ⑤ Now rotate the Zoom Control back to highest magnification and check the focus. Repeat steps ② to ④ to achieve the best focus accuracy.
- ⑥ Look through the left Eyepiece with the left eye and adjust focus using the Left Diopter Adj. Ring similar to steps ① to ⑤.

NSK Niigata seiki Co., Ltd.

5-3-14, Tsukanome, Sanjo, Niigata, Japan, 955-0055
Tel.: +81-256-33-5522 Fax: +81-256-33-5518
MAIL: intl.sales@niigataseiki.co.jp
URL: <http://www.niigataseiki.co.jp>